

Technical Information

iTHERM TM401

Resistance thermometer for hygienic and aseptic applications

Metric version with basic technology
for all standard applications, permanent insert



Applications

- Specially designed for use in hygienic and aseptic applications in the Food & Beverages and Life Sciences industries
- Measuring range: -50 to $+200$ °C (-58 to $+392$ °F)
- Pressure range up to 50 bar (725 psi)
- Protection class: up to IP69K
- Can be used in non-hazardous areas

Head transmitter

All Endress+Hauser transmitters are available with enhanced accuracy and reliability compared to directly wired sensors. Outputs and communication protocol:

Analog output 4 to 20 mA, HART

Your benefits

- Best price-performance ratio and fast delivery time
- User-friendly and reliable from product selection to maintenance
- International certification: hygiene standards as per 3-A, EHEDG, ASME BPE, FDA, TSE Certificate of Suitability
- Wide range of process connections

Table of contents

Function and system design	3	Certificates and approvals	23
iTHERM Hygiene line	3	Hygiene standard	24
Measuring principle	3	Materials in contact with food/product (FCM)	24
Measuring system	4	CRN approval	24
Input	6	Surface cleanliness	24
Measured variable	6	Material resistance	24
Measuring range	6	Ordering information	24
Output	6	Accessories	24
Output signal	6	Device-specific accessories	25
Family of temperature transmitters	6	Communication-specific accessories	26
Power supply	6	Service-specific accessories	26
Wiring diagram for RTD	6	System components	27
Cable entries	8	Supplementary documentation	28
Device plug	8	Brief Operating Instructions (KA)	28
Overvoltage protection	8	Operating Instructions (BA)	28
Performance characteristics	9	Safety Instructions (XA)	28
Reference conditions	9	Functional Safety Manual (FY/SD)	28
Accuracy	9	Registered trademarks	28
Influence of ambient temperature	9		
Self heating	9		
Response time	10		
Calibration	10		
Insulation resistance	10		
Mounting	10		
Orientation	10		
Installation instructions	10		
Environment	14		
Ambient temperature range	14		
Storage temperature	14		
Humidity	14		
Climate class	14		
Degree of protection	14		
Shock and vibration resistance	14		
Electromagnetic compatibility (EMC)	14		
Process	14		
Process temperature range	14		
Thermal shock	14		
Process pressure range	14		
Medium - state of aggregation	15		
Mechanical construction	16		
Design, dimensions	16		
Weight	16		
Material	16		
Surface roughness	17		
Terminal heads	17		
Process connections	19		

Function and system design

iTHERM Hygiene line

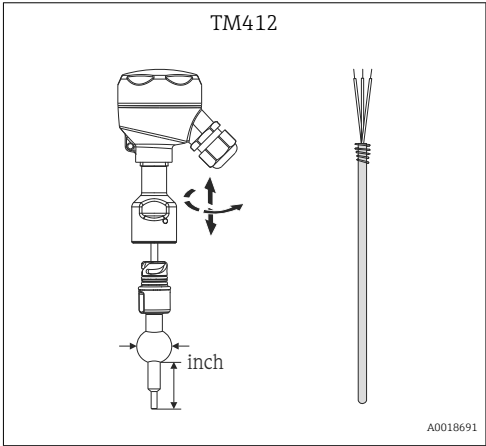
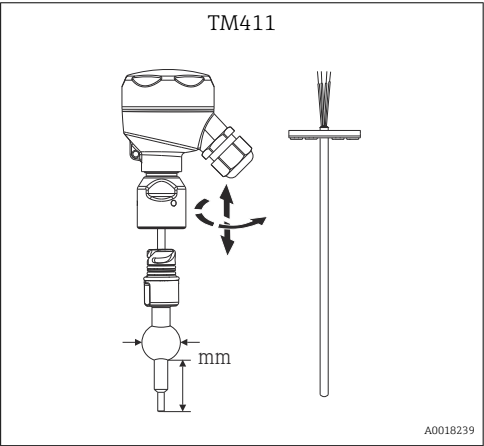
This thermometer is part of the product line of modular thermometers for hygienic and aseptic applications.

Differentiating factors when selecting a suitable thermometer

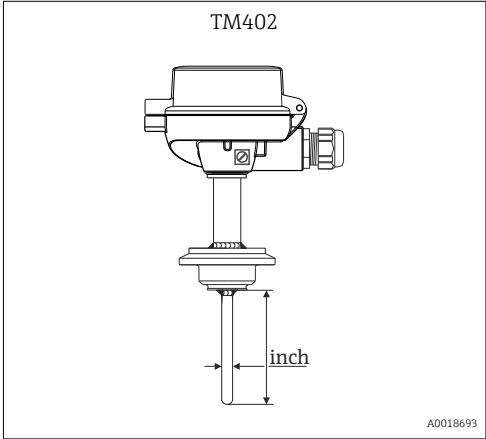
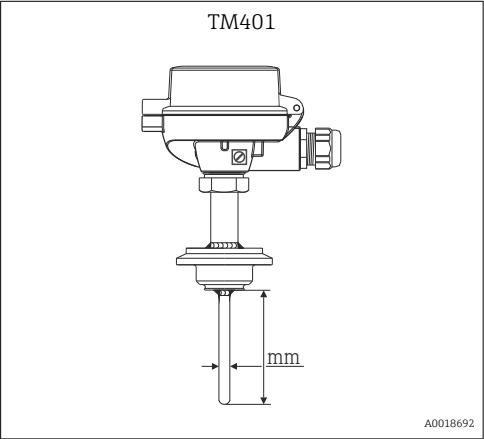
TM4x1	TM4x2
Metric version	Imperial version



TM41x characterizes the device that uses cutting-edge technology, with features such as a replaceable insert, quick-fastening extension neck (iTHERM QuickNeck), vibration-resistant and fast-response sensor technology (iTHERM StrongSens and QuickSens) and approval for use in hazardous areas



TM40x characterizes the device that uses basic technology, with features such as a fixed, non-replaceable insert, application in non-hazardous areas, standard extension neck, low-cost unit



Measuring principle

Resistance thermometer (RTD)

These resistance thermometers use a Pt100 temperature sensor according to IEC 60751. The temperature sensor is a temperature-sensitive platinum resistor with a resistance of 100 Ω at 0 °C (32 °F) and a temperature coefficient $\alpha = 0.003851 \text{ } ^\circ\text{C}^{-1}$.

There are generally two different kinds of platinum resistance thermometers:

- **Wire wound (WW):** Here, a double coil of fine, high-purity platinum wire is located in a ceramic support. This is then sealed top and bottom with a ceramic protective layer. Such resistance thermometers not only facilitate very reproducible measurements but also offer good long-term stability of the resistance/temperature characteristic within temperature ranges up to 600 °C (1 112 °F). This type of sensor is relatively large in size and it is comparatively sensitive to vibrations.
- **Thin film platinum resistance thermometers (TF):** A very thin, ultrapure platinum layer, approx. 1 µm thick, is vaporized in a vacuum on a ceramic substrate and then structured photolithographically. The platinum conductor paths formed in this way create the measuring resistance. Additional covering and passivation layers are applied and reliably protect the thin platinum layer from contamination and oxidation, even at high temperatures.

The primary advantages of thin film temperature sensors over wire wound versions are their smaller sizes and better vibration resistance. A relatively low principle-based deviation of the resistance/temperature characteristic from the standard characteristic of IEC 60751 can frequently be observed among TF sensors at high temperatures. As a result, the tight limit values of tolerance category A as per IEC 60751 can only be observed with TF sensors at temperatures up to approx. 300 °C (572 °F).

Thermocouples (TC)

Thermocouples are comparatively simple, robust temperature sensors which use the Seebeck effect for temperature measurement: if two electrical conductors made of different materials are connected at a point, a weak electrical voltage can be measured between the two open conductor ends if the conductors are subjected to a thermal gradient. This voltage is called thermoelectric voltage or electromotive force (emf.). Its magnitude depends on the type of conducting materials and the temperature difference between the "measuring point" (the junction of the two conductors) and the "cold junction" (the open conductor ends). Accordingly, thermocouples primarily only measure differences in temperature. The absolute temperature at the measuring point can be determined from these if the associated temperature at the cold junction is known or is measured separately and compensated for. The material combinations and associated thermoelectric voltage/temperature characteristics of the most common types of thermocouple are standardized in the IEC 60584 and ASTM E230/ANSI MC96.1 standards.

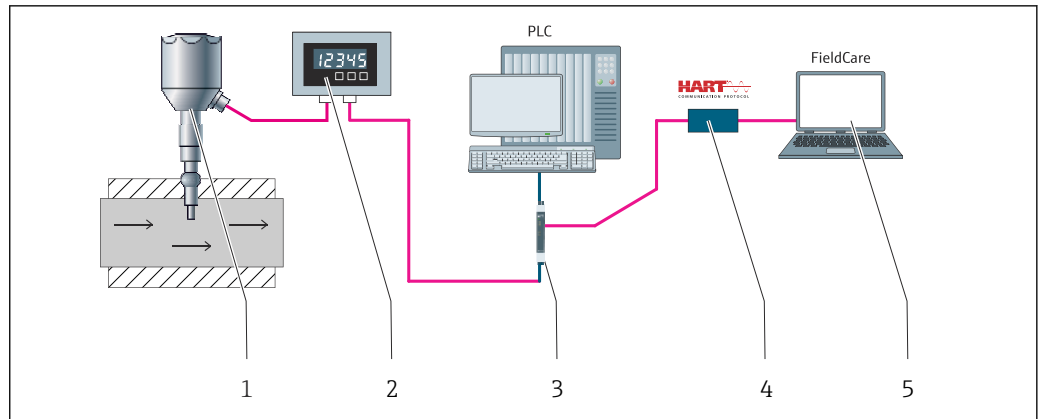
Measuring system

Endress+Hauser offers a complete portfolio of optimized components for the temperature measuring point – everything you need for the seamless integration of the measuring point into the overall facility. This includes:

- Power supply unit/barrier
- Display units
- Overvoltage protection



For more information, see the brochure 'System Components - Solutions for a Complete Measuring Point' (FA00016K/EN)



A0047137

1 Example of application, measuring point layout with additional Endress+Hauser components

- 1 Installed iTHERM compact thermometer with HART communication protocol
- 2 2-wire RIA15 process indicator – The process indicator is looped into the current loop and displays the measuring signal or the HART process variables in digital form. The process indicator does not require an external power supply, as it gets its energy directly from the current loop. More information on this can be found in the Technical Information, under "Documentation".
- 3 Active barrier RN22 – 1- or 2-channel active barrier for the separation of 0/4 to 20 mA standard signal circuits, optionally available as a signal doubler, 24 V DC. HART-transparent. More information on this can be found in the Technical Information, under "Documentation".
- 4 Commubox FXA195 for intrinsically safe HART communication with FieldCare via the USB port.
- 5 FieldCare is a FDT-based plant asset management tool from Endress+Hauser, more details see section 'accessories'. The acquired self-calibration data is stored in the device (1) and can be read using FieldCare. This also enables an auditable calibration certificate to be created and printed.

Input

Measured variable	Temperature (temperature-linear transmission behavior)	
Measuring range	Sensor type	Measuring range
	Pt100 thin-film	-50 to +200 °C (-58 to +392 °F)

Output

Output signal	Generally, the measured value can be transmitted in one of two ways: ▪ Directly-wired sensors - sensor measured values forwarded without a transmitter. ▪ Via all common protocols by selecting an appropriate Endress+Hauser iTEMP temperature transmitter. All the transmitters listed below are mounted directly in the terminal head and wired with the sensory mechanism.
----------------------	--

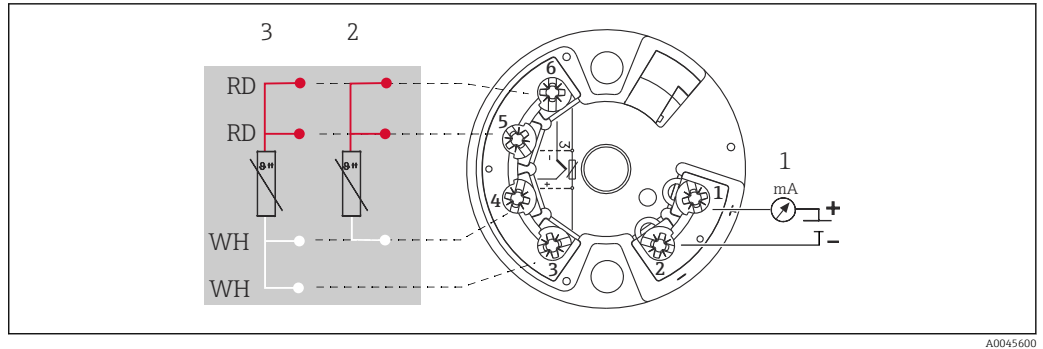
Family of temperature transmitters	Thermometers fitted with iTEMP transmitters are an installation-ready complete solution to improve temperature measurement by significantly increasing accuracy and reliability, when compared to direct wired sensors, as well as reducing both wiring and maintenance costs. PC programmable head transmitters They offer a high degree of flexibility, thereby supporting universal application with low inventory storage. The iTEMP transmitters can be configured quickly and easily at a PC. Endress+Hauser offers free configuration software which can be downloaded from the Endress+Hauser Website. More information can be found in the Technical Information. HART programmable head transmitters The transmitter is a 2-wire device with one or two measuring inputs and one analog output. The device not only transfers converted signals from resistance thermometers and thermocouples, it also transfers resistance and voltage signals using HART communication. It can be installed as an intrinsically safe apparatus in Zone 1 hazardous areas and is used for instrumentation in the terminal head (flat face) according to DIN EN 50446. Quick and easy operation, visualization and maintenance using universal configuration software like FieldCare, DeviceCare or FieldCommunicator 375/475. For more information, see the Technical Information. Advantages of the iTEMP transmitters: ▪ Dual or single sensor input (optionally for certain transmitters) ▪ Unsurpassed reliability, accuracy and long-term stability in critical processes ▪ Mathematical functions ▪ Monitoring of the thermometer drift, sensor backup functionality, sensor diagnostic functions ▪ Sensor-transmitter-matching for dual-channel transmitters, based on Callendar/Van Dusen coefficients
---	--

Power supply



- According to the 3-A Sanitary Standard and the EHEDG, electrical connecting cables must be smooth, corrosion-resistant and easy to clean.
- Grounding or shield connections are possible via special ground terminals on the terminal head.

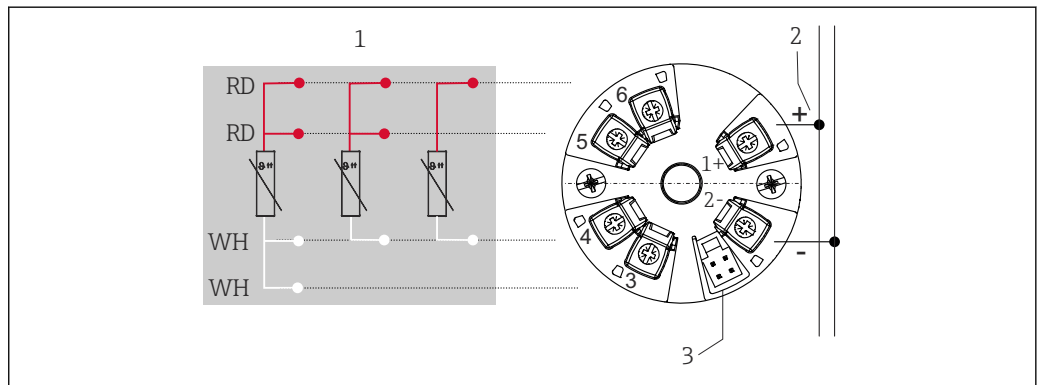
Wiring diagram for RTD	Type of sensor connection
-------------------------------	---------------------------



A0045600

2 Head mounted transmitter TMT18x (single input)

- 1 Power supply for head transmitter and analog output 4 to 20 mA or fieldbus connection
- 2 RTD, 3-wire
- 3 RTD, 4-wire

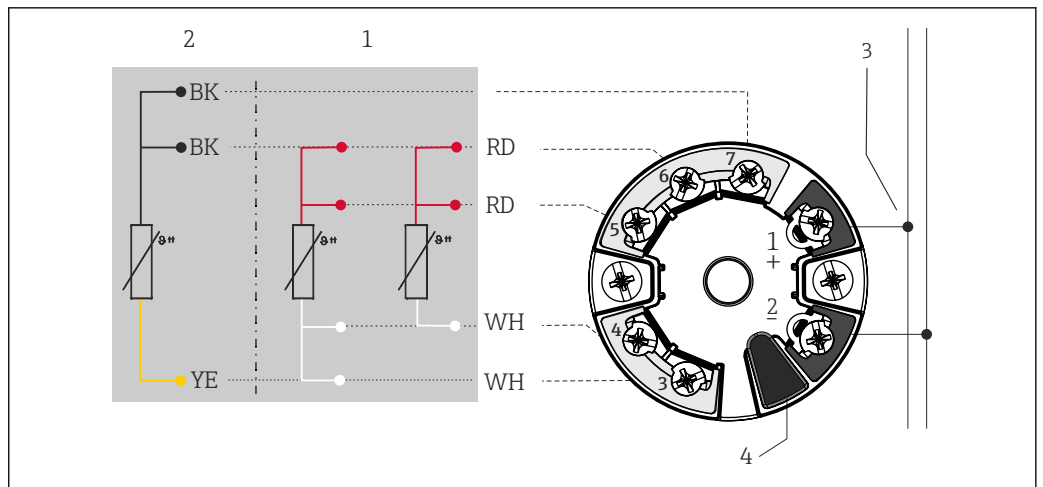


A0047173

3 Head mounted transmitter TMTx1 (single input)

- 1 Sensor input, RTD and Ω : 4-, 3- and 2-wire
- 2 Power supply
- 3 CDI interface, display connection (depends on the head transmitter)

Fitted with push-in terminals if screw terminals are not explicitly selected or a double sensor is installed.

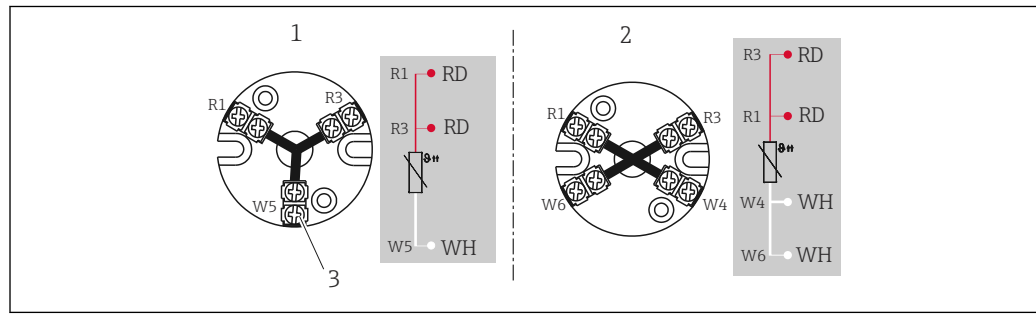


A0045466

4 Head mounted transmitter TMT8x (dual input)

- 1 Sensor input 1, RTD: 4- and 3-wire
- 2 Sensor input 2, RTD: 3-wire
- 3 Power supply or fieldbus connection
- 4 Display connection

Fitted with push-in terminals if screw terminals are not explicitly selected or a double sensor is installed.



A0047086

5 Terminal block mounted

- 1 3-wire, single
2 4-wire, single
3 Outside screw

Cable entries

See "Terminal heads" section

Device plug

PIN assignment of the M12 plugs, connection combinations

Plug	M12 plug, 4-pin			
PIN number	1	2	3	4
Electrical connection (terminal head)				
Flying leads	Not connected (not insulated)			
3-wire terminal block (1x Pt100)	RD	RD	WH	
4-wire terminal block (1x Pt100)			WH	WH
1x TMT 4 to 20 mA or HART	+	i	-	i
PIN position and color code	1 BN 2 GNYE 3 BU 4 GY			

A0018929

Abbreviations

i	RD	WH	BN	GNYE	BU	GY
Insulated ¹⁾	Red	White	Brown	Green-yellow	Blue	Gray

- 1) Wires marked 'i' are not connected and are insulated with heat shrink tubes.

Overvoltage protection

To protect against overvoltage in the power supply and signal/communication cables for the thermometer electronics, Endress+Hauser offers the HAW562 surge arrester for DIN rail mounting and the HAW569 for field housing installation.

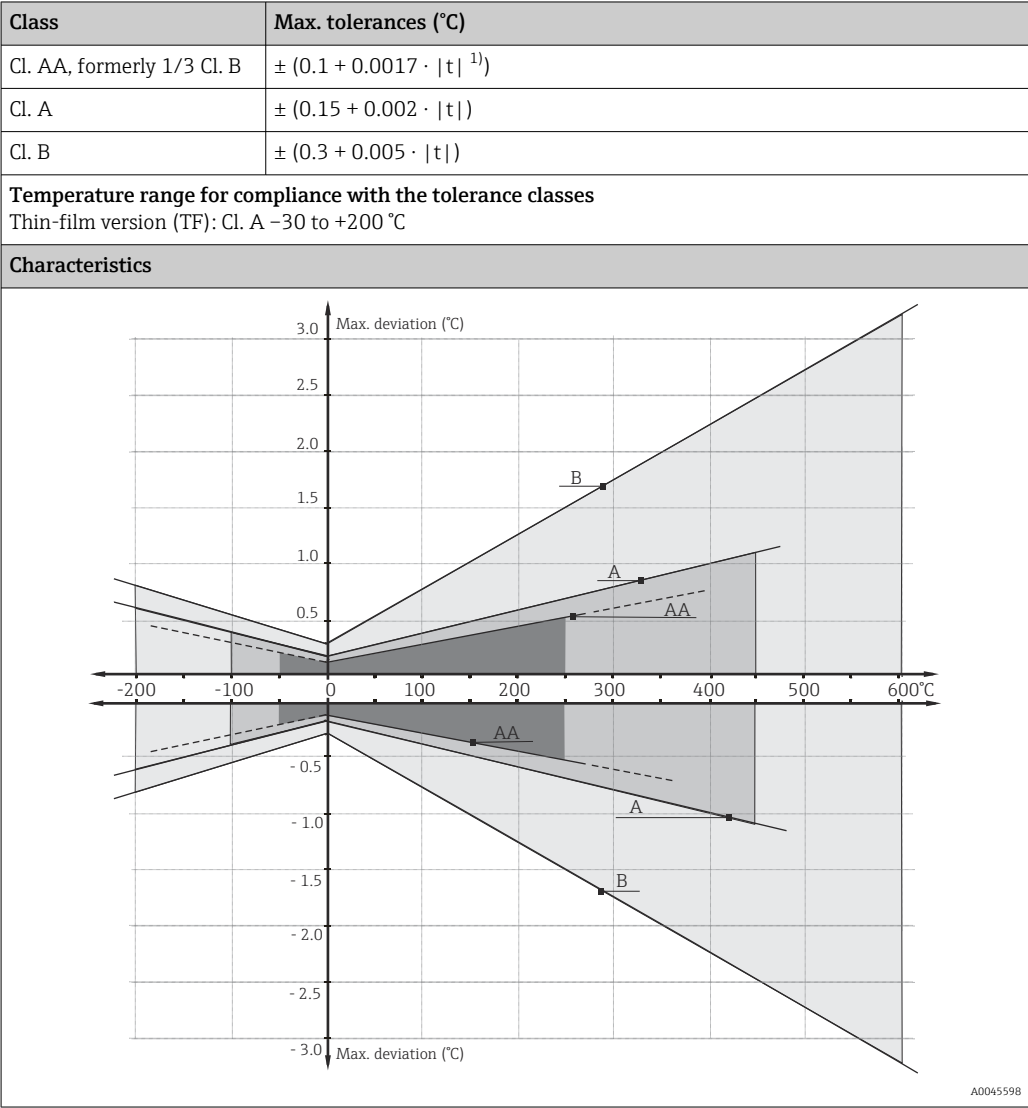


For more information see the Technical Information 'HAW562 Surge arrester' TI01012K and 'HAW569 Surge arrester' TI01013K.

Performance characteristics

Reference conditions These data are relevant for determining the accuracy of the temperature transmitters used. More information on this can be found in the Technical Information of the iTEMP temperature transmitters.

Accuracy RTD resistance thermometer as per IEC 60751



Influence of ambient temperature Depends on the head transmitter used. For details, see the Technical Information.

Self heating RTD elements are passive resistors that are measured using an external current. This measurement current causes a self-heating effect in the RTD element itself which in turn creates an additional measurement error. In addition to the measurement current, the size of the measurement error is also affected by the temperature conductivity and flow velocity of the process. This self-heating error is negligible when an Endress+Hauser iTEMP temperature transmitter (very low measured current) is used.

Response time

Tests have been performed in water at 0.4 m/s (according to IEC 60751) and with a 10 K temperature change.

Pipe diameter	Shape of tip	1x Pt100 thin-film sensor	
		Response time	
		t ₅₀	t ₉₀
Ø6 mm (¼ in)	Straight	5 s	11 s
	Reduced 4.5 mm (0.18 in) x 18 mm (0.71 in)	3.5 s	9 s
Ø8 mm (0.31 in)	Reduced 5.3 mm (0.21 in) x 20 mm (0.79 in)	5 s	10.5 s



Response time without transmitter.

Calibration**Calibration of thermometers**

Calibration involves comparing the measured values of a device under test (DUT) with those of a more precise calibration standard using a defined and reproducible measurement method. The aim is to determine the deviation of the DUT's measured values from the true value of the measured variable. Two different methods are used for thermometers:

- Calibration at fixed-point temperatures, e.g. at the freezing point of water at 0 °C,
- Calibration compared against a precise reference thermometer.

The thermometer to be calibrated must display the fixed point temperature or the temperature of the reference thermometer as accurately as possible. Temperature-controlled calibration baths with very homogeneous thermal values, or special calibration furnaces are typically used for thermometer calibrations. The measuring uncertainty may increase due to heat conduction errors and short immersion lengths. The existing measuring uncertainty is recorded on the individual certificate of calibration. For accredited calibrations in accordance with ISO17025, a measuring uncertainty that is twice as high as the accredited measuring uncertainty is not permitted. If this limit is exceeded, only a factory calibration is possible.

For the device, Endress+Hauser offers standard calibrations at a reference temperature of –50 to +200 °C (–58 to +392 °F) based on the ITS90 (International Temperature Scale). Calibrations in other temperature ranges are available from your Endress+Hauser sales center on request. Calibrations are traceable to national and international standards. The calibration certificate is referenced to the serial number of the thermometer.

Insulation resistance

Insulation resistance ≥ 100 MΩ at ambient temperature, measured between the terminals and the outer jacket with a minimum voltage of 100 V_{DC}.

Mounting

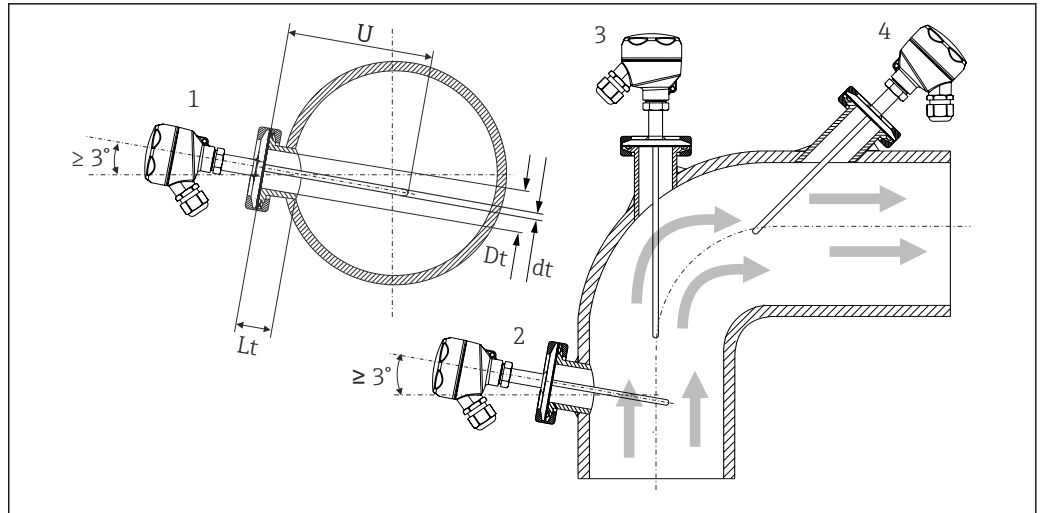
Orientation

No restrictions. However, self-draining in the process must be guaranteed. If there is an opening to detect leaks at the process connection, this opening must be at the lowest possible point.

Installation instructions

The immersion length of the thermometer can influence the accuracy. If the immersion length is too small then errors in the measurement are caused by heat conduction via the process connection and the container wall. If installing in a pipe, the immersion length should ideally correspond to half of the pipe diameter.

- Installation possibilities: pipes, tanks or other plant components
- To minimize the heat conduction error, a minimum immersion length, which corresponds to the calibration, is recommended depending on the type of sensor used.



A0008946

6 Installation examples

- 1, 2 Perpendicular to the flow direction, installed at a min. angle of 3° to ensure self-draining
- 3 On elbows
- 4 Inclined installation in pipes with a small nominal diameter
- U Immersion length

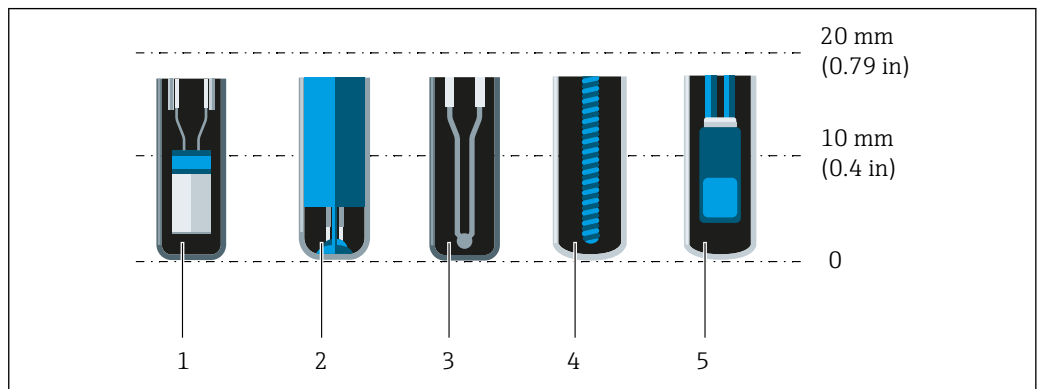
i In the case of pipes with a small nominal diameter, it is advisable for the tip of the thermometer to project well into the process so that it extends past the pipe axis. Installation at an angle (4) could be another solution. When determining the immersion length or installation depth, all the parameters of the thermometer and of the medium to be measured must be taken into account (e.g. flow velocity, process pressure).

i The requirements of the EHEDG and the 3-A Sanitary Standard must be adhered to.

Installation instructions EHEDG/cleanability: $L_t \leq (D_t - d_t)$

Installation instructions 3-A/cleanability: $L_t \leq 2(D_t - d_t)$

Pay attention to the exact position of the sensor element in the thermometer tip.



A0041814

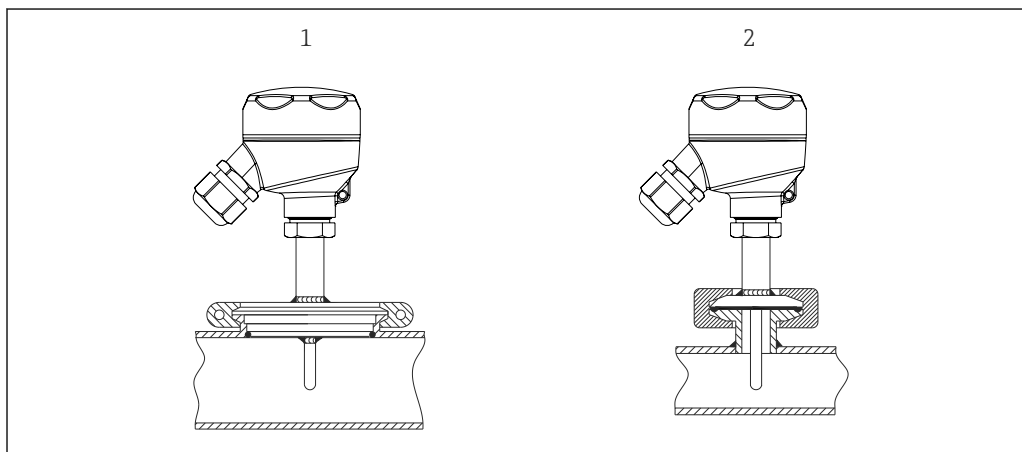
- 1 StrongSens or TrustSens at 5 to 7 mm (0.2 to 0.28 in)
- 2 QuickSens at 0.5 to 1.5 mm (0.02 to 0.06 in)
- 3 Thermocouple (not grounded) at 3 to 5 mm (0.12 to 0.2 in)
- 4 Wire wound sensor at 5 to 20 mm (0.2 to 0.79 in)
- 5 Standard thin-film sensor at 5 to 10 mm (0.2 to 0.39 in)

To keep the influence of heat dissipation to a minimum and to achieve the best possible measurement results, 20 to 25 mm (0.79 to 0.98 in) should be in contact with the medium in addition to the actual sensor element.

This results in the following recommended minimum immersion lengths

- TrustSens or StrongSens 30 mm (1.18 in)
- QuickSens 25 mm (0.98 in)
- Wire wound sensor 45 mm (1.77 in)
- Standard thin-film sensor 35 mm (1.38 in)

It is particularly important to take this into consideration for T-pieces, as the immersion length is very short on account of their design, and the measured error is higher as a result. It is therefore recommended to use elbow pieces with QuickSens sensors.

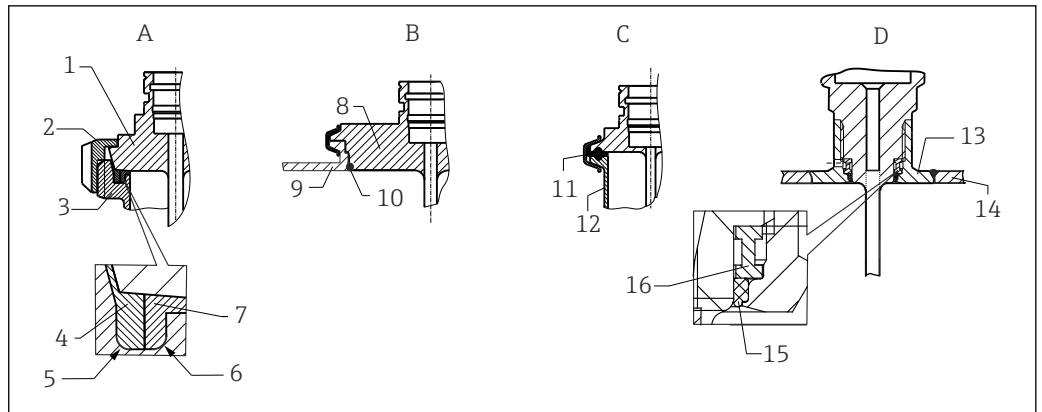


A0018881

 7 Process connections for thermometer installation in pipes with small nominal diameters

1 Varivent® - process connection $D = 50 \text{ mm}$ for DN25 pipes

2 Clamp or micro-clamp



A0040345

8 Detailed installation instructions for hygiene-compliant installation (depends on the version ordered)

A Milk pipe connection according to DIN 11851, only in connection with EHEDG certified and self-centering sealing ring

1 Sensor with milk pipe connection

2 Groove slip-on nut

3 Counterpart connection

4 Centering ring

5 R0.4

6 R0.4

7 Sealing ring

B Varivent® process connection for VARINLINE® housing

8 Sensor with Varivent connection

9 Counterpart connection

10 O-ring

C Clamp according to ISO 2852

11 Gasket seal

12 Counterpart connection

D Process connection Liquiphant-M G1", horizontal installation

13 Weld-in adapter

14 Vessel wall

15 O-ring

16 Thrust collar

NOTICE

The following actions must be taken if a sealing ring (O-ring) or seal fails:

- ▶ The thermometer must be removed.
- ▶ The thread and the O-ring joint/sealing surface must be cleaned.
- ▶ The sealing ring or seal must be replaced.
- ▶ CIP must be performed after installation.

i The counterpieces for the process connections and the seals or sealing rings are not included in the scope of delivery for the thermometer. Liquiphant M weld-in adapters with related seal kits are available as accessories. → 24.

In the case of weld-in connections, exercise the necessary degree of care when performing the welding work on the process side:

1. Use suitable welding material.
 2. Flush-weld or weld with welding radius $\geq 3.2 \text{ mm}$ (0.13 in).
 3. Avoid crevices, folds or gaps.
 4. Ensure the surface is honed and mechanically polished, $R_a \leq 0.76 \text{ }\mu\text{m}$ (30 μin).
1. As a general rule, the thermometers should be installed in such a way that does not impact their ability to be cleaned (the requirements of the 3-A Sanitary Standard must be observed).

2. The Varivent® and Liquiphant-M weld-in adapter and Ingold (+ weld-in adapter) connections enable flush-mounted installation.



For the requirements for installation according to the EHEDG and 3-A Sanitary Standard, see the Operating Instructions for the modular hygienic thermometers.

Operating Instructions BA02023T

Environment

Ambient temperature range

Terminal head	Temperature in °C (°F)
Without mounted head transmitter	Depends on the terminal head used and the cable gland or fieldbus connector, see 'Terminal heads' section
With mounted head transmitter	–40 to 85 °C (–40 to 185 °F)
With mounted head transmitter and display	–20 to 70 °C (–4 to 158 °F)

Terminal head	Temperature in °C (°F)
Without mounted head transmitter	Depends on the terminal head used and the cable gland or fieldbus connector, see 'Terminal heads' section
With mounted head transmitter	–40 to 85 °C (–40 to 185 °F)

Storage temperature

For information, see the ambient temperature.

Humidity

Depends on the transmitter used. If using Endress+Hauser iTEMP head transmitters:

- Condensation permitted as per IEC 60 068-2-33
- Max. rel. humidity: 95% as per IEC 60068-2-30

Climate class

As per EN 60654-1, Class C

Degree of protection

Max. IP69K, depending on the design (terminal head, connector, etc.)

Shock and vibration resistance

The Endress+Hauser inserts meet the requirements of IEC 60751 which specify shock and vibration resistance of 3g in the range from 10 to 500 Hz. The vibration resistance at the measuring point depends on the sensor type and design, see the following table:

Version	Vibration resistance for the sensor tip
Pt100 (TF)	30 m/s ² (3g)

Electromagnetic compatibility (EMC)

Depends on the head transmitter used. For details see the Technical Information.

Process

Process temperature range

Maximum –50 to +200 °C (–58 to +392 °F)

Thermal shock

Thermal shock resistance in CIP/SIP process (temperature increase within 2 seconds from +5 to +130 °C (+41 to +266 °F)).

Process pressure range

The maximum possible process pressure depends on various influencing factors, such as the thermometer design, process connection and process temperature. For information on the maximum

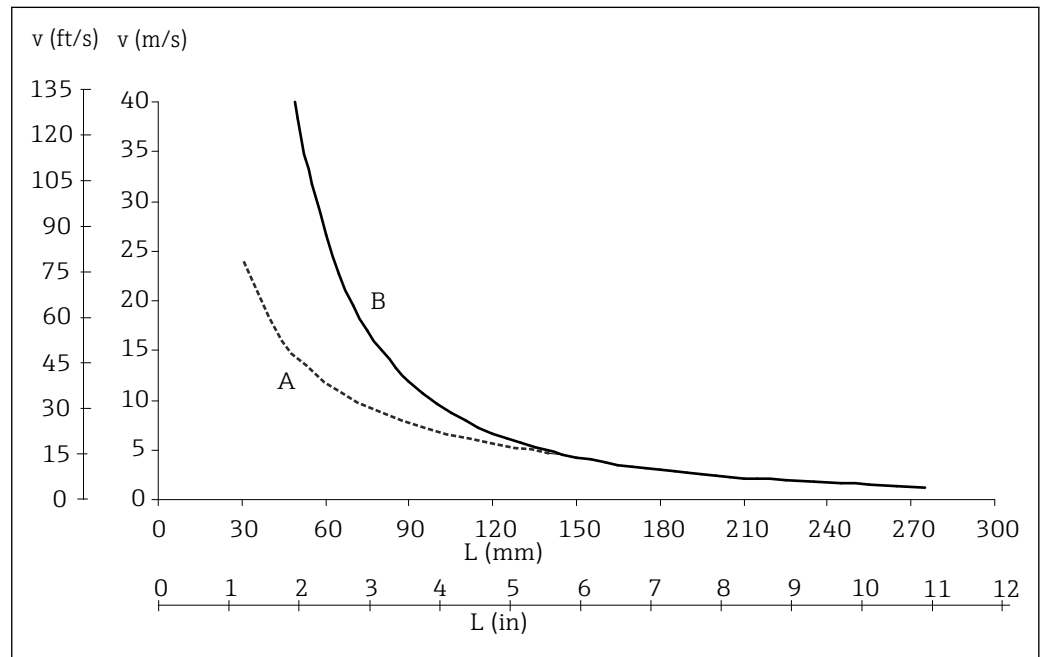
possible process pressures for the individual process connections, see the 'Process connection' section.
→ 19



It is possible to verify the mechanical loading capacity as a function of the installation and process conditions using the online TW Sizing Module for thermowells in the Endress+Hauser Applicator software. See also the 'Accessories' section.

Example of the permitted flow rate as a function of the immersion length and medium

The highest flow velocity tolerated by the thermometer diminishes with increasing thermometer immersion length exposed to the stream of the fluid. The flow velocity is also dependent on the diameter of the thermometer tip, the type of medium being measured, the process temperature and the process pressure. The following figures exemplify the maximum permitted flow velocities in water and superheated steam at a process pressure of 40 bar (580 PSI).



A0008065

9 Permitted flow velocities, thermowell diameter 6 mm (¼ in)

A Medium water at $T = 50\text{ °C}$ (122 °F)

B Medium superheated steam at $T = 400\text{ °C}$ (752 °F)

L Immersion length exposed to flow

v Flow velocity

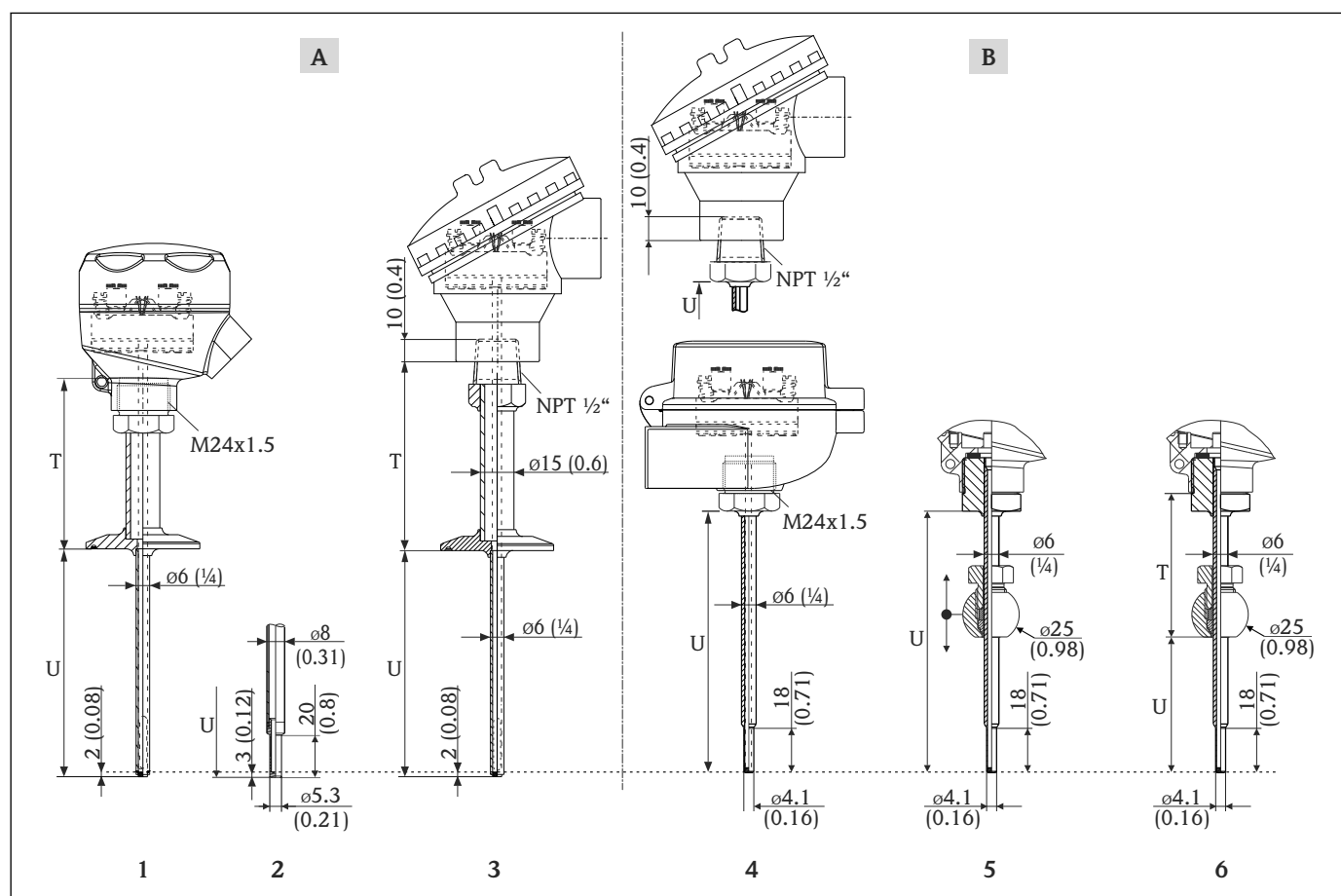
Medium - state of aggregation

Gaseous or liquid (also with high viscosity, e.g. yogurt).

Mechanical construction

Design, dimensions

All dimensions in mm (in).



A0018909

A Version with process connection

B Version without process connection or optionally with compression fitting

1 Thermometer with process connection and M24x1.5 thread to terminal head - shape of tip ø6 mm (0.25 in) straight or

2 Optional shape of tip: ø6 mm (0.25 in) reduced to 5.3 mm (0.21 in) 5.3 mm (0.21 in)

3 Thermometer with process connection and NPT 1/2" thread to the terminal head

4 Thermometer without process connection with M24x1.5 thread (optional NPT 1/2" thread) to terminal head - shape of tip ø6 mm (0.25 in) reduced

5 Thermometer with spherical, movable TK40 compression fitting for weld-in - shape of tip ø6 mm (0.25 in) reduced

6 Thermometer with spherical, fixed TK40 compression fitting for weld-in - shape of tip ø6 mm (0.25 in) reduced

T Length of extension neck (T = 0, for version without process connection or for version with movable compression fitting)

U Immersion length

Weight

0.5 to 2.5 kg (1 to 5.5 lbs) for standard options.

Material

The temperatures for continuous operation specified in the following table are only intended as reference values for use of the various materials in air and without any significant compressive load.

The maximum operating temperatures can be reduced considerably in cases where abnormal conditions such as high mechanical load occur or in aggressive media.

Designation	Short form	Recommended max. temperature for continuous use in air	Properties
AISI 316L (corresponds to 1.4404 or 1.4435)	X2CrNiMo17-13-2, X2CrNiMo18-14-3	650 °C (1 202 °F) ¹⁾	■ Austenitic, stainless steel ■ High corrosion resistance in general ■ Particularly high corrosion resistance in chlorine-based and acidic, non-oxidizing atmospheres through the addition of molybdenum (e.g. phosphoric and sulfuric acids, acetic and tartaric acids with a low concentration) ■ Increased resistance to intergranular corrosion and pitting ■ The wetted part in a protective tube is made of 316L or 1.4435+316L passivated with 3% sulfuric acid.

- 1) Can be used to a limited extent up to 800 °C (1472 °F) for low compressive loads and in non-corrosive media. Contact your Endress+Hauser sales team for further information.

Surface roughness

Values for wetted surfaces:

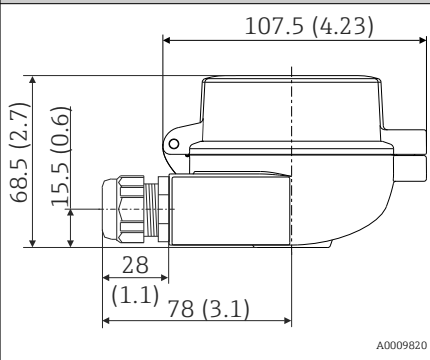
Standard surface, mechanically polished ¹⁾	$R_a \leq 0.76 \mu\text{m}$ (30 μin)
Mechanically polished, buffed ²⁾	$R_a \leq 0.38 \mu\text{m}$ (15 μin)

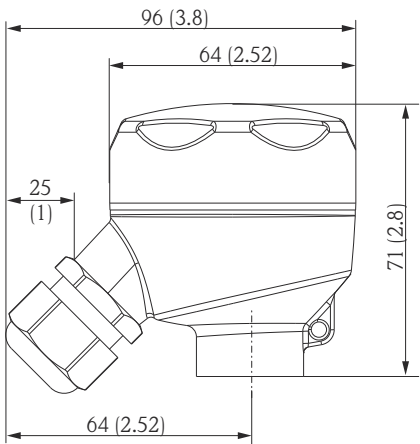
- 1) Or any other finishing method that meets the R_a max
 2) Not compliant with ASME BPE

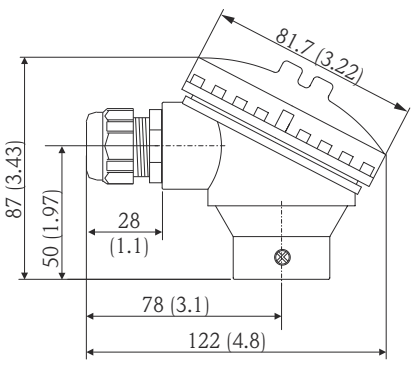
Terminal heads

All terminal heads have an internal shape and size in accordance with DIN EN 50446, flat face and a thermometer connection with a M24x1.5 or ½" NPT thread. All dimensions in mm (in). The sample cable glands in the diagrams correspond to M20x1.5 connections with non-Ex polyamide cable glands. Specifications without head transmitter installed. For ambient temperatures with head transmitter installed, see the 'Environment' section. → 14

As a special feature, Endress+Hauser offers terminal heads with optimized terminal accessibility for easy installation and maintenance.

TA30A	Specification
	■ Degree of protection: ■ IP66/68 (NEMA Type 4x encl.) ■ For ATEX: IP66/67 ■ Temperature: -50 to +150 °C (-58 to +302 °F) without cable gland ■ Material: aluminum, polyester powder coated ■ Seals: silicone ■ Cable entry thread: G ½", ½" NPT and M20x1.5; ■ Protective fitting connection: M24x1.5 ■ Head color: blue, RAL 5012 ■ Cap color: gray, RAL 7035 ■ Weight: 330 g (11.64 oz) ■ Ground terminal, internal and external ■ Available with sensors with 3-A® symbol

TA30R	Specification
 A0018914	■ Degree of protection - standard version: IP69K (NEMA Type 4x encl.) ■ Temperature: -50 to +130 °C (-58 to +266 °F) without cable gland ■ Material: stainless steel 316L, abrasive-blasted or hand-polished ■ Seals: EPDM ■ Cable entry thread ½" NPT and M20x1.5 ■ Weight: 360 g (12.7 oz) ■ Protection armature connection: M24x1.5 or ½" NPT ■ Ground terminal: internal in standard version ■ Not allowed for Class II and III applications ■ Available with 3-A marked sensors

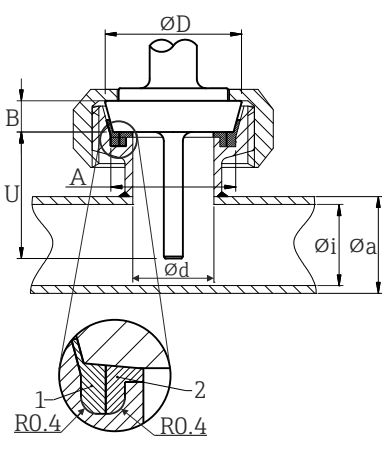
TA30S	Specification
 A0017146	■ Degree of protection: IP65 (NEMA Type 4x encl.) ■ Temperature: -40 to +85 °C (-40 to +185 °F) without cable gland ■ Material: polypropylene (PP), FDA-compliant, seals: O-ring EPDM ■ Cable entry thread: ¾" NPT (with adapter for ½" NPT), M20x1.5 ■ Protective assembly connection: ½" NPT ■ Color: white ■ Weight: approx. 100 g (3.5 oz) ■ Ground terminal: only internal via auxiliary terminal ■ Not allowed for Class II and III applications ■ Available with 3-A marked sensors

Cable glands and fieldbus connectors

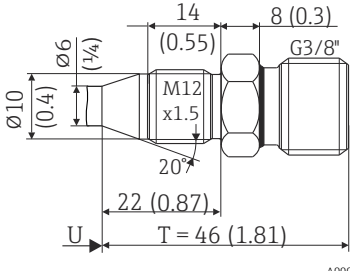
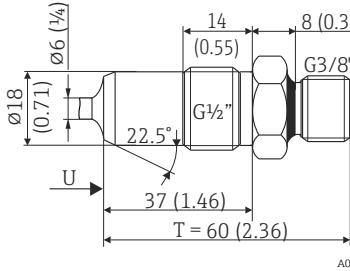
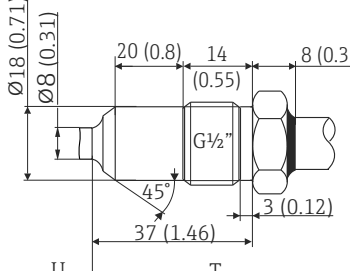
Type	Suitable for cable entry	Degree of protection	Temperature range
Cable gland, polyamide	½" NPT, ¾" NPT, M20x1.5	IP68	-40 to +100 °C (-40 to +212 °F)
	½" NPT, M20x1.5	IP69K	-20 to +95 °C (-4 to +203 °F)
Fieldbus connector (M12, 4-pin)	½" NPT, M20x1.5	IP67, NEMA Type 6	-30 to +90 °C (-22 to +194 °F)

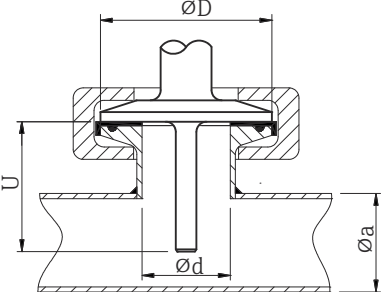
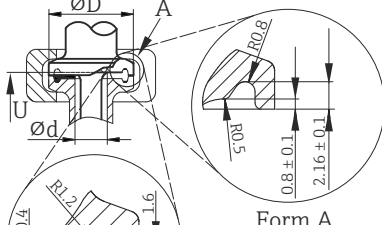
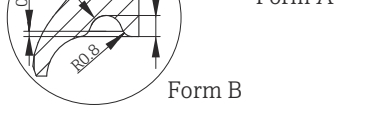
Process connections All dimensions in mm (in).

Releasable process connection

Type						Technical properties
Sanitary connection according to DIN 11851  1 Centering ring 2 Sealing ring A0009561						■ 3-A marked and EHEDG certified (only with EHEDG-certified and self-centering sealing ring). ■ ASME BPE compliance
Version ¹⁾	Dimensions					P _{max.}
	ØD	A	B	Øi	Øa	
DN25	44 mm (1.73 in)	30 mm (1.18 in)	10 mm (0.39 in)	26 mm (1.02 in)	29 mm (1.14 in)	40 bar (580 psi)
DN32	50 mm (1.97 in)	36 mm (1.42 in)	10 mm (0.39 in)	32 mm (1.26 in)	35 mm (1.38 in)	40 bar (580 psi)
DN40	56 mm (2.2 in)	42 mm (1.65 in)	10 mm (0.39 in)	38 mm (1.5 in)	41 mm (1.61 in)	40 bar (580 psi)
DN50	68 mm (2.68 in)	54 mm (2.13 in)	11 mm (0.43 in)	50 mm (1.97 in)	53 mm (2.1 in)	25 bar (363 psi)

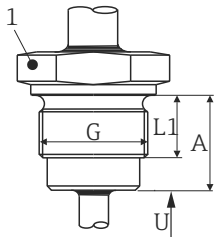
1) Pipes in accordance with DIN 11850

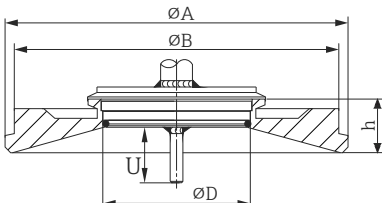
Type	Version	Technical properties
Metal sealing system		
M12x1.5 	G$\frac{1}{2}$" 	Protection pipe diameter 6 mm ($\frac{1}{4}$ in) $P_{max.} = 16$ bar (232 psi)  Maximum torque = 10 Nm (7.38 lbf ft)
-		Protection pipe diameter 8 mm (0.31 in) $P_{max.} = 16$ bar (232 psi)  Maximum torque = 10 Nm (7.38 lbf ft)

Type	Version	Dimensions		Technical properties	Conformity
	ϕd^1	ϕD	ϕa		
Clamp according to ISO 2852    Form A: In compliance with ASME BPE Type A Form B: In compliance with ASME BPE Type B and ISO 2852 Reference: A0009566	Microclamp ²⁾ DN8-18 (0.5"-0.75") ³⁾ , Form A	25 mm (0.98 in)	-	$P_{max.} = 16$ bar (232 psi), depends on clamp ring and suitable seal 3-A marked	ASME BPE Type A
	Tri-clamp DN8-18 (0.5"-0.75"), Form B		-		Based on ISO 2852 ⁴⁾
	Clamp DN12-21.3, Form B	34 mm (1.34 in)	16 to 25.3 mm (0.63 to 0.99 in)		ISO 2852

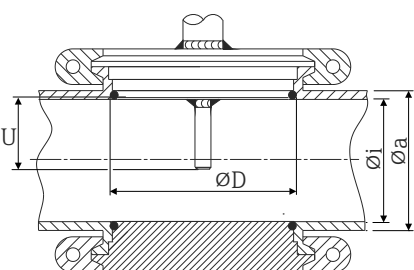
Type	Version	Dimensions		Technical properties	Conformity
	ϕd ¹⁾	ϕD	ϕa		
	Clamp DN25-38 (1"-1.5"), Form B	50.5 mm (1.99 in)	29 to 42.4 mm (1.14 to 1.67 in)	■ $P_{max.} = 16$ bar (232 psi), depends on clamp ring and suitable seal ■ 3-A marked and EHEDG certified (in connection with Combifit seal) ■ Can be used with "Novaseptic Connect (NA Connect)" which enables flush-mount installation	ASME BPE Type B; ISO 2852
	Clamp DN40-51 (2"), Form B	64 mm (2.52 in)	44.8 to 55.8 mm (1.76 to 2.2 in)		ASME BPE Type B; ISO 2852

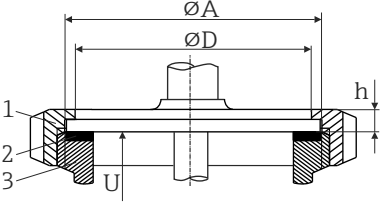
- 1) Pipes in accordance with ISO 2037 and BS 4825 Part 1
2) Microclamp (not in ISO 2852); no standard pipes
3) DN8 (0.5") only possible with protection pipe diameter = 6 mm (¼ in)
4) Groove diameter = 20 mm

Type	Version G	Dimensions			Technical properties
		L1 thread length	A	1 (SW/AF)	
Thread according to ISO 228 (for Liquiphant weld-in adapter) 	G¾" for FTL20/31/33 adapter	16 mm (0.63 in)	25.5 mm (1 in)	32	■ $P_{max.} = 25$ bar (362 psi) at max. 150 °C (302 °F) ■ $P_{max.} = 40$ bar (580 psi) at max. 100 °C (212 °F) ■ Information about hygienic compliance in connection with FTL31/33/50 adapter see TI00426F
	G¾" for FTL50 adapter				
	G1" for FTL50 adapter	18.6 mm (0.73 in)	29.5 mm (1.16 in)	41	

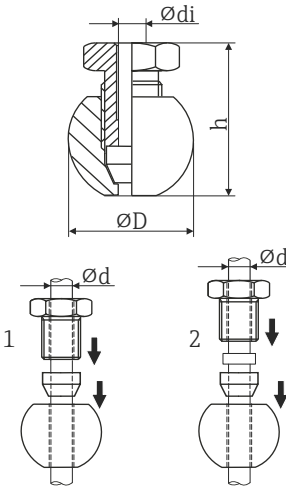
Type	Version	Dimensions				Technical properties	
		ϕD	ϕA	ϕB	h	$P_{max.}$	
Varivent® 	Type F	50 mm (1.97 in)	145 mm (5.71 in)	135 mm (5.31 in)	24 mm (0.95 in)	10 bar (145 psi)	■ 3-A marked and EHEDG certified ■ ASME BPE compliance
	Type N	68 mm (2.67 in)	165 mm (6.5 in)	155 mm (6.1 in)	24.5 mm (0.96 in)		

i The VARINLINE® housing connection flange is suitable for welding into the conical or torispherical head in tanks or containers with a small diameter (≤ 1.6 m (5.25 ft)) and up to a wall thickness of 8 mm (0.31 in).

Type				Technical properties	
Varivent® for VARINLINE® housing for installation in pipes  A0009564				■ 3-A marked and EHEDG certified ■ ASME BPE compliance	
Version	Dimensions			P _{max.}	
	ØD	Øi	Øa		
Type N, according to DIN 11866, series A	68 mm (2.67 in)	DN40: 38 mm (1.5 in)	DN40: 41 mm (1.61 in)	DN40 to DN65: 16 bar (232 psi)	
		DN50: 50 mm (1.97 in)	DN50: 53 mm (2.1 in)		
		DN65: 66 mm (2.6 in)	DN65: 70 mm (2.76 in)		
		DN80: 81 mm (3.2 in)	DN80: 85 mm (3.35 in)	DN80 to DN150: 10 bar (145 psi)	
		DN100: 100 mm (3.94 in)	DN100: 104 mm (4.1 in)		
		DN125: 125 mm (4.92 in)	DN125: 129 mm (5.08 in)		
		DN150: 150 mm (5.9 in)	DN150: 154 mm (6.06 in)		
Type N, according to EN ISO 1127, series B	68 mm (2.67 in)	38.4 mm (1.51 in)	42.4 mm (1.67 in)	42.4 mm (1.67 in) to 60.3 mm (2.37 in): 16 bar (232 psi)	
		44.3 mm (1.75 in)	48.3 mm (1.9 in)		
		56.3 mm (2.22 in)	60.3 mm (2.37 in)		
		72.1 mm (2.84 in)	76.1 mm (3 in)	76.1 mm (3 in) to 114.3 mm (4.5 in): 10 bar (145 psi)	
		82.9 mm (3.26 in)	42.4 mm (3.5 in)		
		108.3 mm (4.26 in)	114.3 mm (4.5 in)		
Type N, according to DIN 11866, series C	68 mm (2.67 in)	OD 1½": 34.9 mm (1.37 in)	OD 1½": 38.1 mm (1.5 in)	OD 1½" to OD 2½": 16 bar (232 psi)	
		OD 2": 47.2 mm (1.86 in)	OD 2": 50.8 mm (2 in)		
		OD 2½": 60.2 mm (2.37 in)	OD 2½": 63.5 mm (2.5 in)		
Type N, according to DIN 11866, series C	68 mm (2.67 in)	OD 3": 73 mm (2.87 in)	OD 3": 76.2 mm (3 in)	OD 3" to OD 4": 10 bar (145 psi)	
		OD 4": 97.6 mm (3.84 in)	OD 4": 101.6 mm (4 in)		

Type	Version	Dimensions			Technical properties
		ϕD	ϕA	h	
SMS 1147  A0009568 1 Coupling nut 2 Sealing ring 3 Counterpart connection	DN25	32 mm (1.26 in)	35.5 mm (1.4 in)	7 mm (0.28 in)	P _{max.} = 6 bar (87 psi)
	DN38	48 mm (1.89 in)	55 mm (2.17 in)	8 mm (0.31 in)	
	DN51	60 mm (2.36 in)	65 mm (2.56 in)	9 mm (0.35 in)	
 The counterpart connection must fit the sealing ring and fix it in place.					

Without process connection (for compression fitting)

Type	Version	Dimensions			Technical properties ¹⁾
		Ødi	ØD	h	
Compression fitting TK40 for weld-in  1 Movable 2 Fixed A0018912	Spherical Material of sealing taper PEEK or 316L Thread G $\frac{1}{4}$ "	6.3 mm (0.25 in) for protection pipe diameter Ød = 6 mm (0.236 in)	25 mm (0.98 in)	33 mm (1.3 in)	■ P_{max.} = 10 bar (145 psi), T_{max.} = +150 °C (+302 °F) for PEEK material, tightening torque = 10 Nm ■ P_{max.} = 50 bar (725 psi), T_{max.} = +200 °C (+392 °F) for 316L material, tightening torque = 25 Nm ■ PEEK compression fitting is EHEDG tested, 3-A marked

1) All the pressure specifications apply for cyclic temperature load



The 316L compression fittings can only be used once due to deformation. This applies to all the components of the compression fittings! A replacement compression fitting must be attached in a different position (grooves in the protection pipe). PEEK compression fittings must never be used at a temperature that is lower than the temperature present when the compression fitting is secured. This is because the fitting would no longer be leak-tight as a result of heat contraction of the PEEK material.

SWAGELOCK or similar fittings are strongly recommended for higher requirements.

Minimum extension neck lengths, dependent on respective process connection

Process connection	Extension next length T
■ None ■ Compression fitting, movable	Predefined (cannot be selected, T = 0)
■ Thread according to ISO 228 ■ Compression fitting, fixed ■ Metal sealing system	≥82 mm (3.23 in)
■ Clamp according to ISO 2852 ■ Sanitary connection according to DIN 11851 ■ Varivent[®] ■ SMS 1147	≥55 mm (2.17 in)

Certificates and approvals

Current certificates and approvals that are available for the product can be selected via the Product Configurator at www.endress.com:

1. Select the product using the filters and search field.
2. Open the product page.
3. Select **Configuration**.

Hygiene standard	■ EHEDG certification, type EL CLASS I. EHEDG certified/tested process connections. →  19 ■ 3-A authorization no. 1144, 3-A Sanitary Standard 74-07. Listed process connections. →  19 ■ ASME BPE, certificate of conformity can be ordered for indicated options ■ FDA-compliant ■ All surfaces in contact with the medium are free of animal derived ingredients (ADI/TSE) and do not contain any materials derived from bovine or animal sources.
Materials in contact with food/product (FCM)	The materials of the thermometer in contact with food/product (FCM) comply with the following European regulations: ■ (EC) No. 1935/2004, Article 3, paragraph 1, Articles 5 and 17 on materials and articles intended to come into contact with food. ■ (EC) No. 2023/2006 on good manufacturing practice for materials and articles intended to come into contact with food. ■ (EU) No. 10/2011 on plastic materials and articles intended to come into contact with food.
CRN approval	The CRN approval is only available for certain thermowell versions. These versions are identified and displayed accordingly during the configuration of the device. Detailed ordering information is available for your nearest sales organization www.addresses.endress.com or in the Download Area under www.endress.com : 1. Select the country 2. Select Downloads 3. In the search area: select Approvals/approval type 4. Enter the product code or device 5. Start the search
Surface cleanliness	Free from oil and grease for O ₂ applications, optional
Material resistance	Material resistance (including housing) to the following cleaning agents/disinfectants from the company Ecolab: P3-topax 66, P3-topactive 200, P3-topactive 500 and P3-topactive OKTO as well as demineralized water.

Ordering information

Detailed ordering information is available for your nearest sales organization www.addresses.endress.com or in the Product Configurator under www.endress.com :

1. Click Corporate
2. Select the country
3. Click Products
4. Select the product using the filters and search field
5. Open the product page

The Configuration button to the right of the product image opens the Product Configurator.



Product Configurator - the tool for individual product configuration

- Up-to-the-minute configuration data
- Depending on the device: Direct input of measuring point-specific information such as measuring range or operating language
- Automatic verification of exclusion criteria
- Automatic creation of the order code and its breakdown in PDF or Excel output format
- Ability to order directly in the Endress+Hauser Online Shop

Accessories

Various accessories, which can be ordered with the device or subsequently from Endress+Hauser, are available for the device. Detailed information on the order code in question is available from your

local Endress+Hauser sales center or on the product page of the Endress+Hauser website:
www.endress.com.

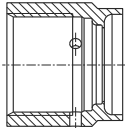
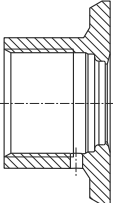
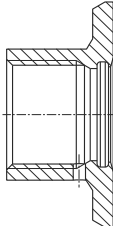
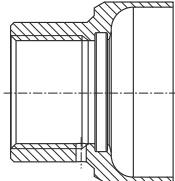
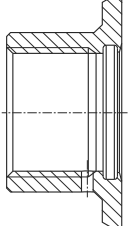
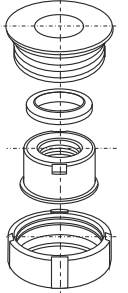
Device-specific accessories

Accessories	Description
Welding boss with sealing taper (metal - metal) A0006621 A0018236	Welding boss for G$\frac{1}{2}$" and M12x1 thread Metal-sealing; conical Material of wetted parts: 316L/1.4435 Max. process pressure 16 bar (232 PSI) Order number: ■ 71424800 (G$\frac{1}{2}$") ■ 71405560 (M12x1)
Dummy plug A0045726 1 Size across flats SW22	Dummy plug for G$\frac{1}{2}$" or M12x1 conical metal-sealing welding boss Material: SS 316L/1.4435 Order number: ■ 60022519 (G$\frac{1}{2}$") ■ 60021194 (M12x1)
Weld-in adapter for Ingold process connection (OD25 mm (0.98 in)x46 mm (1.81 in) A0008956	Material of wetted parts: 316L/1.4435 Weight: 0.32 kg (0.7 lb) Adapter for Ingold process connection with 3.1 material certificate, order number: 71531585 Adapter for Ingold process connection, order number: 71531588 O-ring seal set ■ Silicone O-ring in accordance with FDA CFR 21 ■ Maximum temperature: 230 °C (446 °F) ■ Order number: 60018911

Weld-in adapter



For more information about order codes and hygienic compliance of the adapters and spare parts, see Technical Information (TI00426F).

Weld-in adapter	 A0008246	 A0008251	 A0008256	 A0011924	 A0008248	 A0008253
	G 3/4", d=29 for pipe-mounting	G 3/4", d=50 for vessel-mounting	G 3/4", d=55 with flange	G 1", d=53 without flange	G 1", d=60 with flange	G 1" adjustable
Material	316L (1.4435)	316L (1.4435)	316L (1.4435)	316L (1.4435)	316L (1.4435)	316L (1.4435)
Roughness μm (μin) process side	≤ 1.5 (59.1)	≤ 0.8 (31.5)	≤ 0.8 (31.5)	≤ 0.8 (31.5)	≤ 0.8 (31.5)	≤ 0.8 (31.5)

i Maximum process pressure for the weld-in adapters:

- 25 bar (362 PSI) at maximum 150 °C (302 °F)
- 40 bar (580 PSI) at maximum 100 °C (212 °F)

Communication-specific accessories

Configuration kit TXU10	Configuration kit for PC-programmable transmitter with setup software and interface cable for PC with USB port Order code: TXU10-xx
Commubox FXA195 HART	For intrinsically safe HART communication with FieldCare via the USB port.  For details, see "Technical Information" TI00404F
Wireless HART adapter SWA70	Is used for the wireless connection of field devices. The WirelessHART adapter can be easily integrated into field devices and existing infrastructures, offers data protection and transmission safety and can be operated in parallel with other wireless networks with minimum cabling complexity.  For details, see Operating Instructions BA061S
Fieldgate FXA320	Gateway for the remote monitoring of connected 4-20 mA measuring devices via a Web browser.  For details, see "Technical Information" TI00025S and Operating Instructions BA00053S

Service-specific accessories

Accessories	Description
Applicator	Software for selecting and sizing Endress+Hauser measuring devices: ■ Calculation of all the necessary data for identifying the optimum measuring device: e.g. pressure loss, accuracy or process connections. ■ Graphic illustration of the calculation results Administration, documentation and access to all project-related data and parameters over the entire life cycle of a project. Applicator is available: Via the Internet: https://portal.endress.com/webapp/applicator

Accessories	Description
Configurator	Product Configurator - the tool for individual product configuration ■ Up-to-the-minute configuration data ■ Depending on the device: Direct input of measuring point-specific information such as measuring range or operating language ■ Automatic verification of exclusion criteria ■ Automatic creation of the order code and its breakdown in PDF or Excel output format ■ Ability to order directly in the Endress+Hauser Online Shop The Configurator is available on the Endress+Hauser website at: www.endress.com -> Click "Corporate" -> Select your country -> Click "Products" -> Select the product using the filters and search field -> Open product page -> The "Configure" button to the right of the product image opens the Product Configurator.
DeviceCare SFE100	Configuration tool for devices via fieldbus protocols and Endress+Hauser service protocols. DeviceCare is the tool developed by Endress+Hauser for the configuration of Endress+Hauser devices. All smart devices in a plant can be configured via a point-to-point or point-to-bus connection. The user-friendly menus enable transparent and intuitive access to the field devices.  For details, see Operating Instructions BA00027S
FieldCare SFE500	FDT-based plant asset management tool from Endress+Hauser. It can configure all smart field units in your system and helps you manage them. By using the status information, it is also a simple but effective way of checking their status and condition.  For details, see Operating Instructions BA00027S and BA00065S
Accessories	Description
W@M	Life cycle management for your plant W@M offers assistance with a wide range of software applications over the entire process: from planning and procurement to the installation, commissioning and operation of the measuring devices. All the relevant information is available for every measuring device over the entire life cycle, such as the device status, device-specific documentation, spare parts etc. The application already contains the data of your Endress+Hauser device. Endress+Hauser also takes care of maintaining and updating the data records. W@M is available: Via the Internet: www.endress.com/lifecyclemanagement

System components

Accessories	Description
RIA15 field indicator	The process indicator is looped into the current loop and displays the measuring signal or the HART process variables in digital form. The process indicator does not require an external power supply, as it gets its energy directly from the current loop.  For details, see "Technical Information" TI01043K
RN22	1- or 2-channel active barrier for separation of 0/4 to 20 mA standard signal circuits, optionally available as a signal doubler, 24 V DC. HART-transparent.  For details, see "Technical Information" TI01515K
RNS221	Supply unit for powering two 2-wire measuring devices solely in the non-Ex area. Bidirectional communication is possible via the HART communication jacks.  For details, see "Technical Information" TI00081R and Brief Operating Instructions KA00110R

Supplementary documentation

The following document types are available in the Downloads section of the Endress+Hauser website (www.endress.com/downloads):



For an overview of the scope of the associated Technical Documentation, refer to the following:

- *W@M Device Viewer* (www.endress.com/deviceviewer): Enter the serial number from the nameplate
- *Endress+Hauser Operations App*: Enter the serial number from the nameplate or scan the matrix code on the nameplate

Brief Operating Instructions (KA)

Guide that takes you quickly to the 1st measured value

The Brief Operating Instructions contain all the essential information from incoming acceptance to initial commissioning.

Operating Instructions (BA)

Your reference guide

These Operating Instructions contain all the information that is required in various phases of the life cycle of the device: from product identification, incoming acceptance and storage, to mounting, connection, operation and commissioning through to troubleshooting, maintenance and disposal.

Safety Instructions (XA)

Depending on the approval, the following Safety Instructions (XA) are supplied with the device. They are an integral part of the Operating Instructions.



The nameplate indicates the Safety Instructions (XA) that are relevant to the device.

Functional Safety Manual (FY/SD)

Depending on the SIL approval, the Functional Safety Manual (FY/SD) is an integral part of the Operating Instructions and applies in addition to the Operating Instructions, Technical Information and ATEX Safety Instructions.



The different requirements that apply for the protective function are described in the Functional Safety Manual (FY / SD).

Registered trademarks

HART®

Registered trademark of the FieldComm Group, Austin, Texas, USA



71562674

www.addresses.endress.com
